

4.2. Changing Chosen Wavelengths (metallic type only)

1. Select “Laser” and enter.
2. Select the wavelength you wish to use and enter. If you wish to save this wavelength as the startup default, press save before exiting.
3. If you wish to modify one of the wavelengths, select the closest wavelength, press the right navigation key.
4. Now, using the up/down keys to change the numbers and the right/left keys to move to the next number, modify to the wavelength you want.
5. When finished, press “save” and exit. If you only want a temporary change, exit without “save”.

4.3. To Set Type of Laser Being Used

From main measurement screen, select “Laser” and enter. Choose the desired laser type or wavelength. If you want this one to be the startup wavelength, press save before exiting.

4.4. Setting Startup Configuration

1. Turn on the Nova II with the sensor connected, select “Menu” and enter. Select “Configure”.
2. Now, with the navigation keys select “Startup Mode”, select “power” or “energy” as the startup mode. Press enter to exit.
3. Now, with the navigation keys select “Display Mode”. Press the enter key and select the type of screen you wish to start up with. Press save and exit. If you do not want to save this as the startup mode but only change for now, exit without save. (This choice is saved in the instrument and not the sensor so the Nova II will start up from the screen selected for all sensors).
4. Now select “Energy Range”, press enter and choose the manual energy range you wish to be the default.
5. Now select “Average” to select what period you wish to average power over or select the default, “none”.
6. If the sensor is a diffuser sensor, now select diffuser in or out.
7. Now select “Max Pulse Len” and choose the shortest pulse length setting longer than your laser’s pulse length.

Warning: Incorrect readings will result if pulse length is not set up correctly.

8. Select “Laser” and select the laser type you wish to be the default.
9. Press “save” to save all present settings then “exit” to get back to the main measurement screen.

Note: Some sensors have only the long pulse setting for all pulse lengths. In that case N/A appears. For further details on configuration see Sections 6.2 and 6.3 of the manual.

4.5. Energy or Average Power Measurement

Warning:

Do not exceed maximum sensor limits for power, energy, power density and energy density as listed in tables 5 and 6 section 10.2 of the manual, Sensor Specifications. Otherwise, there is a risk of damaging the absorber.

With the pyroelectric sensor, you have been supplied a test slide with the same coating as on your pyroelectric detector. You can also obtain this slide from your dealer. You should use this slide to test the damage threshold with your laser pulses. If the slide is damaged, then either enlarge your beam or lower the laser energy until damage is no longer seen.

4.6. To choose Energy or Average Power Measurement

To switch from the main energy measurement screen to power measurement, press “power”. To switch from the main power measurement screen to energy measurement, press “energy”.



NOVA II

Quick Reference

1 Getting Started

The Nova II is equipped with “soft keys”. That is, the functions of the keys change as indicated by the legend above each key. It is also equipped with 4 navigation keys and a round enter key in the center.

To connect sensor to the Nova II meter:

Insert the 15 pin D type connector of the measuring sensor cable into the socket marked “Head Input” on the rear panel of the Nova II meter.

To switch the Nova II on:

To switch on, briefly press the on/off/backlight switch (leftmost key). The unit will switch on, and the display will appear. Note that the sensor must be plugged in before the unit is switched on. To switch the backlight on or off, press the on/off/backlight switch briefly again. To switch the Nova II off, press the on/off/backlight switch and hold it for ~ 3 seconds until the display blanks.

If you wish to save the current Nova II configuration, use the “configure” function. (See section 2.3.2.).

To set the display modes:

1. Disconnect the sensor and switch off then on again. “Head Disconnected” will appear.
2. Select “Display Mode” and enter. Choose the desired mode, press save. The Nova II will now be in this display mode for all sensors.
3. Select “Date” and press the enter button. Using the navigation keys set to the correct date. Press save.
4. Select “Time” and enter. Using the navigation keys, set to the correct time. Press save.
5. For adjusting RS232 baud rate and An. Output voltage see the main manual.

Nova II Quick Reference



Ophir Optronics Solutions LTD.
Science Based Industry Park P.O.B. 45021, Jerusalem 9145001, Israel
Phone +972-2-5484444
<https://www.ophiropt.com/photonics>
rev 1.45-5/17.Feb.19/be Ophir P/N 1J06036

To zero instrument:

1. Disconnect the sensor, turn off then on again. “Head Disconnected” will be displayed.
2. Make sure instrument is not in an electrically noisy environment and is undisturbed.
3. Press “zero” and “start”. Wait until “zeroing completed” appears. For more details, see main manual section 3.5.5.

2 Thermal Sensors

2.1. Use of Nova II with Thermal Type Sensors

To set type of laser being used:

1. While the Nova II is off, plug in the sensor then switch it on again.
2. Press the navigation keys until “laser” is highlighted. Press the enter key and select the appropriate laser wavelength.
3. If you want to save this new wavelength as the startup default, press “save”.
4. Return to main screen by pressing the enter key.

To choose manual or automatic ranging in power measurement:

1. Select “Range” with the navigation keys. Press the enter key.
2. Select the appropriate manual range or AUTO.
3. If you want to save this new range as the startup default, press “save”.
4. Press the enter key to return to the main measurement screen.

To choose power or energy measurement:

To go from the main power measurement screen to the energy measurement screen, press the lower left soft key “energy”. To go from the energy measurement screen to power measurement, press “power”.

2.2. Setting and Saving the Startup Configuration

1. From the main screen, select “Menu” and enter. Select “Configure”.
2. Now, with the navigation keys select “Startup Mode”, press the enter key and select “power” or “energy” as the startup mode. Press Enter to return to screen.
3. Now, select “Display Mode”. Press enter and select the startup type of screen. Save if you wish or exit by pressing enter (This choice is saved in the instrument so the Nova II will start up from the screen selected for all sensors).
4. Now select “Power Range”, press enter and choose the manual power range you wish to be the default or choose autorange.
5. Now select “Average” to select what period you wish to average power over or select the default, “none”. See section 4.4.2.4 of the main manual for details.
6. Now select “Energy Range” and choose the default energy range.
7. Now select “Threshold” if you wish to change the energy threshold. Select “Laser” to select the laser type you wish to be the default. Press save if you want to save this configuration, then press exit to return to the main measurement screen.
Further details on configuration in section 4.3. of the main manual.

2.3. Power or Single Shot Energy Measurement:

Warning:

Do not exceed maximum sensor limits for power, energy, power density and energy density as listed in tables 5 and 6 in section 10.2. of the main manual. Otherwise, there is a risk of damaging the absorber.

To use the Nova II to measure laser power:

Verify that the display shows “Menu: Power” in the upper right hand side of the display. If the display shows “Menu: Energy” and units of J, mJ etc. then press “power” to switch to the power measurement mode.

To select an analog needle type display for the Nova II:

Select “Menu” and enter. Select “Needle” and enter. The display will now show a simulated analog needle type display. In order to expand the display, press “Expand”.

To expand the bargraph scale $\pm 5x$ about the present reading:

1. From the main power measurement screen press the “Zoom” button.
2. Press the “Zoom” button again to return to the ordinary bargraph.

To subtract background and set current reading to zero:

1. From the main power measurement screen press the “Offset” button.
2. Press “Offset” again to cancel. See Section 4.4.2.2. of the main manual for full details.

To use the Nova II to fine-tune laser power:

1. From the main power measurement screen select “Menu” and enter. Now select “Tune” and enter.
2. Set the percentage range of the power scale to be displayed by repeatedly pressing the $\pm 50\%$ key.
3. Now select the horizontal sweep time you desire by selecting “Graph” and choosing the time scale.
4. You may now use the screen to tune the laser power. See Section 4.4.2.3. of the main manual for full details.
5. Press “power” to exit to the main measurement screen.

To use the Nova II to measure laser energy:

1. Verify that the display shows “energy” in the upper right hand side of the display. If the display shows “power” and units of W, mW etc. then press “energy” to switch to the energy measurement mode.
2. The energy mode is manual ranging. Select “Range”, enter then use the navigation keys to select the proper range. Press save if you want this range to be the startup default. The correct range is the lowest one that is larger than the pulse energy measured. Note that the range will be displayed in the lower right corner of the main display.
3. When the Nova II screen flashes “ready”, on and off, fire the laser. See Section 4.5 of the main manual for full details.

3 Photodiode Sensors

3.1. Selecting Chosen Wavelengths

1. While the Nova II is off, plug in the sensor then switch it on again. From the main measurement screen, press “Laser” to select the correct laser wavelength. If you want to save this new wavelength as the startup default, press “save” before exiting.
2. If the wavelength you want is not among the wavelengths in the six wavelength list and you want to change or add a wavelength, then move to the wavelength you wish to change or add. Press the right navigation key.
3. Using the up/down keys to change each number and the right left keys to move to the next number, key in the wavelength you want. Press enter twice. If you wish to save this new wavelength as a startup default, press “save”, then press enter.
4. Follow the directions on configuration in section 3.2. below.

To choose manual or automatic ranging or dBm in power measurement:

1. Press “Range” and then select the appropriate manual range, autorange or dBm (logarithmic scale).
2. Press enter to return to the main measurement screen. If you wish to save the selected range as the startup default, press “save” before exiting.

3.2. Setting the Startup Configuration

1. Select menu and press enter. Select “configure” and enter.
2. Select “Filter” and enter. Choose filter in or out to be the default.
3. Now select “Display Mode” and choose which screen you want the instrument to start up in. Press enter. (Note that this is an instrument setting and will be valid for all sensors.)
4. Now select “Power Range” and choose the manual power range you wish to be the default or choose autorange or dBm.
5. Now select “Average” and choose which period to average power over or leave the default “NONE”.
6. Now select “Laser” to choose the startup laser wavelength. When you are finished, if you want to save this as default startup, press save. If you only want to change for now, exit without saving. For further details on configuration see Section 5.3 of the main manual.

3.3. Power Measurement

Warning:

Do not exceed maximum sensor limits for power, energy, power density and energy density as listed in tables 5 and 6 section 10.2 of the manual, Sensor Specifications. Otherwise, there is a risk of damaging the absorber.

To expand the bargraph scale $\pm 5x$ about the present reading:

1. From the main power measurement screen, press the “Zoom” button.
2. Press the “Zoom” button again to return to the ordinary bargraph.

To subtract background and set current reading to zero:

1. From the main power measurement screen press the “Offset” button.
2. Press “Offset” again to cancel. See Section 5.4.5. of the main manual for details.

To use the Nova II to fine-tune laser power:

1. From the main power measurement screen select “Menu” and enter. Now select “Tune” and enter.
2. Set the percentage range of the power scale to be displayed by repeatedly pressing the $\pm 50\%$ key.
3. Now select the horizontal sweep time you desire by selecting “Graph” and choosing the time scale.
4. You may now use the screen to tune the laser power. Press “power” to exit to the main measurement screen. See Section 4.4.2.3. of the manual for details.

4 Pyroelectric or Photodiode Energy Sensors

4.1. Zeroing Instrument Against Sensor

For most accurate calibration, you should zero the pyroelectric sensor against the Nova II it is being used with. Proceed as follows: Make sure the sensor is in a quiet environment and not subject to pulsed radiation. Plug sensor into Nova II and turn on. Press “menu”, select “zero” and enter. Press “start”. When “zeroing completed” appears, press “save”.